

Charcoal filter replacement




Attention the exhausted filter is contaminated with harmful substances what formaldehyde or its substitutive.

The hood ecos trim advance needs a periodic substitution of the filters in how much the duration of the filters is influenced by different variable as: the temperature, the damp, the speed of aspiration and the volume of the inhaled vapours.

Diapath guarantees only the complete functionality of the aspiration if filters are installed planned by Diapath s.p.a., in how much you filter different characteristic have they would damage the components climbed on board of the cloak and they would put to risk the health of the operator.

To perform the substitution of the filter to active carbons is owed::


- Abbassare il vetro fino a fine corsa
- Extinguish the tool and therefore disconnect the power supply
- Lift up the door of the upper part
- unscrew the panel that contains the filters

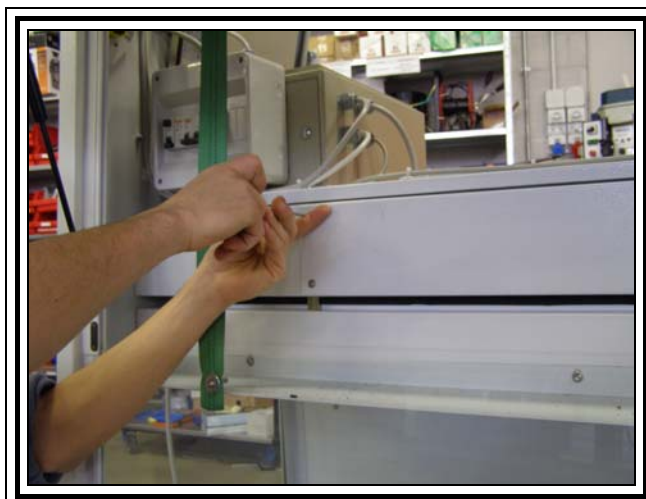

- replace the contaminated filters with those new



- screw the panel that contains the filters
- lower the counter

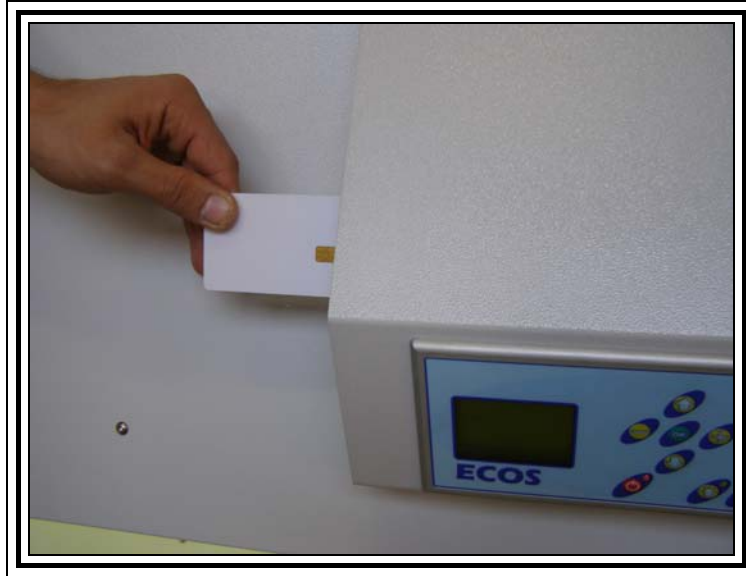

- turn on the cloak and to insert the card Badge in the situated reader in the vain one of the hood